



**TECHNICAL
ADVANTAGES**

Authorised **OHAUS** Australia Dealer
www.techadv.com.au
sales@techadv.com.au

RANGER COUNT 3000

RANGER COUNT 4000

2nd Scale Interface Instruction Manual

INTRODUCTION

This Second Scale interface Kit is for use with the Ohaus product Ranger Count 3000. Through this kit you can connect additional platform with Ranger Count 3000.


KIT CONTENTS

- 2nd Scale PC Board
- Pressure Button
- Screws (4)
- Instruction Manual


INTERFACE INSTALLATION

Install the Interface module on the underside of the balance as shown.


1. Remove the red interface options cover. Break off the plastic blocker from the option cover.

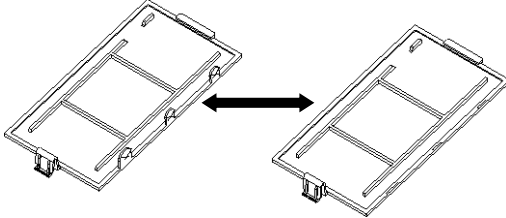

2. Install the Interface PCB. Push it through the pins and secure it with the 4 screws.

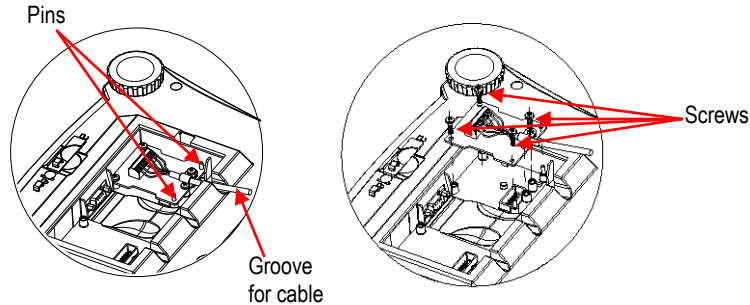


3. Connect a cable.

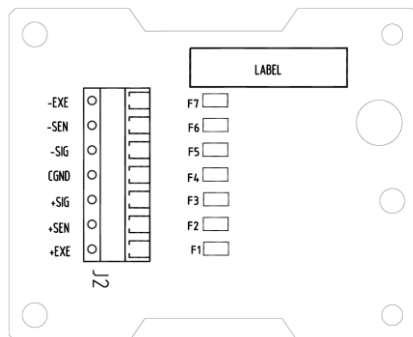

4. Place the cord in the groove and put the red interface option cover back on.

SETUP

After the main platform is connected to the 2nd weighing platform, switch on the Dual Scale (D.Scale) menu item in the SETUP menu (see menu structure below).

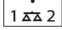Then set the CAP2 and Grad2 in the SETUP menu and do a calibration. Press and hold the  key to switch between the main platform and the secondary platform.

CAL

Zero2
Span1
Span2
Lin1
Lin2
GEO
End

SETUP

Reset
D.Scale
Cap2
Grad2
Pwr. Un
Zero
A.Opt
A.Tare
Bp.Opt
Bp.Sig
Bp.Key
Accum
End

Note: Make sure to perform a calibration of the scale before using the 2nd weighing platform. Make sure the weighing unit is in kg or lb when using high capacity platform. (If display shows error8.6, long press [**Units**] key to switch unit to kg or lb.)

ACCESSORIES

For a complete listing of Ohaus printers and other accessories, contact Ohaus Corporation or visit www.Ohaus.com

COMPLIANCE

This accessory has been tested and complies with the approvals listed in the applicable Instruction Manual.

Disposal



In conformance with the European Directive 2002/96 EC on Waste Electrical and Electronic Equipment (WEEE) this device may not be disposed of in domestic waste. This also applies to countries outside the EU, per their specific requirements.

Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. If you have any questions, please contact the responsible authority or the distributor from which you purchased this device.

Should this device be passed on to other parties (for private or professional use), the content of this regulation must also be related.

For disposal instructions in Europe, refer to www.ohaus.com/weee.



OHAUS Corporation
7 Campus Drive
Parsippany, NJ 07054, USA
Tel: +1 973 377 9000
Fax: +1 973 944 7177

With offices worldwide

www.ohaus.com



* 3 0 0 6 4 4 8 1 *

P/N 30064481B © 2013 Ohaus Corporation, all rights reserved

Printed in China