



www.techadv.com.au sales@techadv.com.au

SF40A PRINTER Instruction Manual



CONTENTS

0	Overview1				
1	Introdu	ction	2		
		Description Features			
2	Startup	Procedure	3		
	2.1	Safety Measures	3		
	2.2	Power Supply	3		
	2.3	Printer setup	4		
		2.3.1 Printer setup using the accompanying software2.3.2 Printer setup using printer only			
	2.4	Installing the Printer	8		
		2.4.1 Connecting the printer to the instrument2.4.2 Switching the printer on2.4.3 Inserting paper	8		
3	Keystro	oke Functions	9		
4	Menu		10		
5	Printer	Test	13		
6	Mainte	nance	13		
	6.1	Changing the Paper Roll	13		
	6.2	Replacing the Ribbon	13		
	6.3	Cleaning	14		
	6.4	Disposal	14		
7	Trouble	shooting	15		
8	Recomr	nended Settings	17		
9	9 Consumables and Spare Part18				
1(10 RS232C Interface18				
1	11 Technical Data19				

Overview

Controls Connections Image: Control of the state of the

5 Interface connector

Control panel

Functions

Key SF40A	
• Zeroing /Taring the • Abort – Exit	e balance
 Printout of date an Scroll down menu 	
Paper feed Scroll up menu op	tion
 Printing of stable v values Opening the menu- select next menu Save settings 	-
• Totaling applicatio	n
Statistics application	on

1 Introduction

1.1 Description

The SF40A is an advanced dot matrix printer for OHAUS instruments. The printer has Statistics and Totalization functions. The SF40A also meets the requirements of modern quality assurance systems such as (GLP, GMP, ISO9001 etc.).

1.2 Features

- Supports all 13 languages present in Ohaus instruments.
- Statistics and Totalization functions.
- Zero / Tare function.
- Real-time clock function.
- The printout meets the requirements of modern quality assurance systems (GLP, GMP, ISO9001, etc.)
- Complies with the requirements of measurement traceability by precisely documenting the origin of data, as well as calibration procedures.
- Meets the requirement of archiving measurement results by providing a thermally stable and light resistant printout.
- Easy setup with the software provided with the printer.

2. Startup Procedure

2.1 Safety Measures

The Printer has been tested for the connections and intended purposes documented in these Operating Instructions. However, it is your responsibility to perform the necessary tests related to the methods and purposes of its intended use. Therefore, you should observe the following safety measures.

- The printer must be used indoors only, and never in hazardous areas.
- The electronics of the printer are protected against the ingress of splashing water and dust. However, since the printer housing is not watertight, the printer should not be used in the presence of liquids.
- Comply with cleaning instruction (see chapter 6.3), protect power plug against wet conditions.
- Never open the printer housing. If you have any problems with your printer, please contact your responsible OHAUS service.
- If the equipment is used in a manner not specified by the manufacturer, the protection provided by the equipment may be impaired.

2.2 Power Supply

Your printer is supplied with an AC adapter with a country-specific power cable. The power supply is suitable for all line voltages in the range: 100 - 240 VAC, 50 - 60 Hz (for exact specifications, see chapter 9).



Attention

- First, check the local line voltage is in the range 100 240 VAC, 50 - 60 Hz and whether the power plug fits your local power supply connection. If this is not the case, on no account connect the printer or the AC adapter to the power supply, but contact the responsible OHAUS dealer.
- Only plug the adapter into a socket which is grounded.

Important

- \triangle
- Before operating, check all cables for damage.
- Guide the cables so that they cannot become damaged or interfere with the weighing process.
- Make sure that the AC adapter cannot come into contact with liquids
- The power plug must always be accessible.

2.3 Printer Setup

The default setting of the printer is as below, if you want change some of them you need set up before use it.

Header:	No Header 🔹	Baud rate:	9600 🔻
Date Format:		Data bits:	8 🗸
Time Format:	24:MM:SS	Parity:	None
Char Set:	CP1252 Window Latin1	Stop bits:	1
Balance Feature:	Balance Feature Enable 🔹 🔻	Flow control:	Xon/Xoff 🔹
O/T Settings:	Perform Zero 🔹		

There are two ways to change the default setting:

- Use the accompanying software
- Operate on the printer only.


2.3.1 Printer setup using accompanying software

- 1) Attach the interconnect cable to the printer interface on the computer and to the printer.
- 2) Connect the printer to the main power supply.
- 3) Use the main switch <I/O> to power on.
- 4) Insert the CD into the computer and install the printer software.
- 5) Run the software and set up the RS232 parameter of computer, then click the "Connect" button.

CO	/ Settings			
Please config the COM port setting.				
	Port:	COM1	Baud rate:	9600 💌
	Data bits:	8 🗸	Parity:	None 🔻
	Stop bits:	1	Flow control:	Xon/Xoff 🔹
		Connect	Disconnect	

SF40A Printer

6) The following window will appear if connection is successful, click OK to continue.

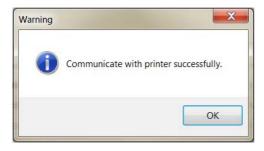

7) The printer Menu setting is now activated and all the settings can be changed.

lenu Setting				
Date Time: 💿 K	eep Printer Time	Current System Time	Custom	2012/08/01 09:51:35
Header:	No Header	•	Baud rate:	9600 🔻
Date Format:	DD.MMM YYYY	•	Data bits:	8
Time Format:	24:MM:SS	•	Parity:	None
Char Set:	CP1252 Window Latin1	•	Stop bits:	1 •
Balance Feature:	Balance Feature Enable	•	Flow control:	Xon/Xoff 🔹
O/T Settings:	Perform Zero	•		
Description: CP1252 languages.	covers Western Europea	n languages such as Eng	lish, German, Frei	nch, Spanish, Italian and other
Factory Reset			Get	Status Save Status

- a. Date Time: Set the date and time for the printer
 - Keep Printer Time: keep the existing date and time in the printer
 - Current System Time: Get the date and time from your computer
 - · Custom: Set the date and time manually

- b. Header: Set the header of the printing content
 - No header: No header for printing content
 - Date/Time: Header with date and time
 - Date/Time/Balance: Header with date, time and instrument ID
- c. Date Format: Set date format
- d. Time Format: Set Time Format
- e. Char Set: Set the Code page for different languages
 - CP 1252 Window Latin1: Covers Western European languages such as English, German, French, Spanish and Italian.
 - CP 1251Window Cyrillic: Covers languages that use the Cyrillic alphabet such as Russian, Bulgarian and Serbian Cyrillic.
 - CP437 MS-DOS Latin US: Covers Western European languages such as English, German, French, Spanish and Italian.
 - CP852 MS-DOS Latin2: Covers Central European languages that use Latin script such as Polish Bosnian, Czech, Hungarian, Croatian, Romanian, Serbian and Slovak.
 - CP860 MS-DOS Portuguese: Covers Portuguese language


f. Balance Feature :

- 1) Enable/Disable a function: 0/T, Statistics, Totalization
- 2) Printer connection status: LED light will either blink or stay lit.
 - Balance Feature Enabled: 0/T, Statistics and Totalization are enabled. The printer connection status LED light will blink if the connection with balance is lost.
 - Balance Feature Disable: 0/T, Statistics, Totalization functions are disabled. The printer connection status LED will be lit even if the connection with balance is lost.

SF40A Printer

- Perform Zero: Perform Zero function when pressing the button.
- Perform Tare: Perform Tare function when pressing the button.
- h. **Baud rate, Data bits, Parity, Stop bits, Flow control:** Set the parameters of RS232 setting.


The settings should be the same as the instrument to be connected. For the instrument RS232 settings, please see the instrument manual.

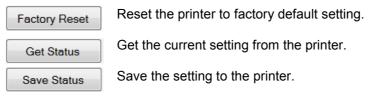

8) Click "Save Status" button to save the setting and click Exit to finish the setup.

2.3.2 Printer setup using printer only

- 1) Connect the printer to the main power supply.
- 2) Use the main switch <I/O> to power on.
- 3) Press and hold the « button to enter the menu list, the printer will print

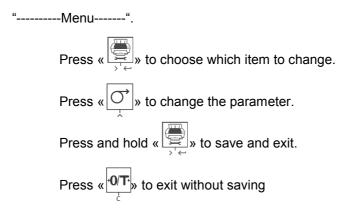

The printer will print the content to show the result of every operation. For total menu structure, please see chapter 4.

2.4 Installing the printer

2.4.1 Connecting the printer to the instrument

Before connecting, make sure the instrument and printer have the same RS232 parameter setting. For instrument Baud rate settings, please see the related manual.

- Attach the interconnect cable to the printer interface of the instrument and to the printer.
- Switch the instrument on.

2.4.2 Switching the printer on




- Connect the printer to the mains power.
- Use the main switch <I/O> to power on.
- RS232 interface only: The pilot lamp will flash until an instrument is connected. If the lamp continues to flash see chapter 7.

2.4.3 Inserting Paper




- Remove paper cover (pull on the back) (1).
- Insert the paper rotary axis through the core of the paper roll (2).
- Feed paper through the slot in the printing unit in the direction of the arrow (3).
- Press and hold until enough paper has been fed (4).
- Replace the paper cover (1).

3 Keystroke Functions

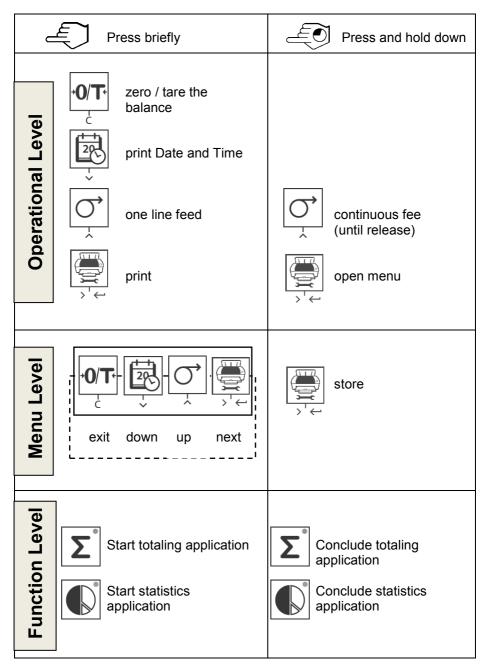

EN-10

4 Menu

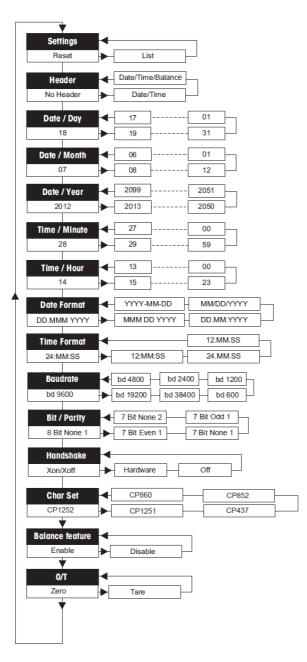

SF40A Printer

Example Printout:

```
-----MENU-----
Settings:
 * List
Header:
 * No Header
Date Day:
   31
Date Month:
   07
Date Year:
   2012
Time Minute:
   0.0
Time Hour:
  18
Date Format:
  * YYYY-MM-DD
Time Format:
  * 24:MM:SS
Baudrate:
  * 9600
Bit/Parity:
  * 8 Bit None 1
Handshake:
  * Xon/Xoff
Char Set:
  * CP1252 Windows Latin 1
Balance Feature:
  * Enable
->0/T<- Setting:
  * Zero
```

* Factory settings

Applications

Press briefly					
Statistics	Statistics Application ¹⁾ Totaling Application ²⁾				
	→ Start "Statistics application" (LED on)	Σ	→ Start "Totaling application" (LED on)		
	→ Add a new sample and press « .».	Σ°	→ Add a new sample and press « ∑ [®] ».		
	→ If needed print subtotal press « → ↔ »	Č,	→ If needed print subtotal press « → · ← »		
	→ Print and end "Statistic application" (LED off).	Σ° C	→ Print and end "Totaling application" (LED off).		

Important notes

Unit change during the application is not permitted.

Maximal 999 sample. ¹⁾ Sample weight must be in the range of 70% - 130% of the current average value.

Minimum dynamic weight must be 100d. ²⁾ Minimum weights + 1 d.

5 Printer Test

The printer is equipped with a self-test that automatically prints out the character set of the printer (standard IBM/DOS).

Starting the test	→	Press the « » key and power on the printer. The printer prints out the current settings.
Terminating the test	→	Power off the printer.

6 Maintenance

Under normal conditions the printer requires practically no maintenance. The OHAUS service department is available if servicing is required. Please ask your responsible OHAUS dealer for details.

6.1 Changing the Paper Roll

Insertion of a new roll of paper is described in chapter 2.3. Please see chapter 9 (Consumables) for paper roll order number. **Note:** You may obtain the paper rotary axis from the old paper roll.

6.2 Replacing the Ribbon



- Remove the paper cover (1)
- Pull paper out of printing unit (2)



- To remove the ribbon cartridge, press in direction of arrow (3).
- Insert the new ribbon, and if necessary, take up slack with tensioning wheel (4). See chapter 9 (Consumables) for the ribbon cartridge order number.
- Insertion of a new roll of paper is described in chapter 2.4.

EN-14

6.3 Cleaning

Since the printer housing is made of top grade, resistant materials, all commercially available mild cleaning agents may be used (see safety precautions in chapter 2.1).

6.4 Disposal



In conformance with the European Directive 2002/96/EC on Waste Electrical and Electronic Equipment (WEEE) this device may not be disposed of in domestic waste. This also applies to countries outside the EU, per their specific requirements.

Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. If you have any questions, please contact the responsible authority or the distributor from which you purchased this device.

Should this device be passed on to other parties (for private or professional use), the content of this regulation must also be related.

Disposal instructions in Europe are available online at europe.ohaus.com/europe/en/home/support/weee.aspx.

Thank you for your contribution to environmental protection.


7 Troubleshooting

Problem / Message	Possible cause	Correction
Printout unreadable	 Ribbon worn or entangled. Service life of printing unit at an end. 	 Change ribbon cartridge or tense the ribbon properly. Please contact OHAUS service.
Pilot lamp does not light	No power.Power supply fault.	 Switch printer on. Please contact OHAUS service.
Pilot lamp flashes	 No connection printer device. Printer did not find the settings of the connected instrument. 	 Make sure that the printer and the instrument is connected. Switch the printer off/on. Alternatively change the settings of the printer or the connected instrument. (see chapter 2.4)
Timeout Over Load	 Data transfer has not completed in the last 45 s. Balance in 	 Move balance to more stable location. Check the balance. Check the sample
	 Verload range. Zeroing out of the zero range. 	 weight. Change the zero range of the balance. Change the setting of the printer from zero to tare.
Under Load	 Balance in underload range. Weighing pan is not in position. 	Check the balance.
Transmission Error	 Instrument does not understand printer instruction. 	 Check interface parameter settings on printer and instrument.

Problem / Message	Possible cause	Correction
Not Executable	Instrument cannot execute printer instruction.	Check the balance.
Date/Time Not Set Memory Error	Printer error messages.	 Check the instrument. Reset the printer. Set date/time Back-up battery depleted Please contact OHAUS service.
Memory full	Statistics or Totaling Memory is full.	Start a new Statistics or Totaling application.
Load/Unload Weight	 No sample or the same sample is on the weighing pan. 	 Put a new sample or remove it and put it again on the pan.
Unit change not allowed!	 Changed unit during Statistics or Totaling application. 	 Reset the balance. Change the unit back to the first one.
Out of range	 Sample weight is out of range 	 Sample weight must be in the range of 70% - 130% of the current average value.


Connected Device	Device Settings	Printer Settings
DV, AV, PA, AR, SP, TA, NV, NVL, NVT, MB35/45, MB23/25, T71P, T32XW, T32WE, T51 XW, T71XW	Printer Mode Baud rate: 9600 Bit / Parity: 8 bit None 1 Handshake: Xon/Xoff	Factory Settings
EX	Printer Mode Baud rate: 9600 Bit / Parity: 8 bit None 1 Handshake: Xon/Xoff	Factory Settings
Russian	Printer Mode Baud rate: 9600 Bit / Parity: 8 bit None 1 Handshake: Xon/Xoff Language: ANSI/WIN Russian	CP1251 Windows Cyrillic
Polish	Printer Mode Baud rate: 9600 Bit / Parity: 8 bit None 1 Handshake: Xon/Xoff Language: ANSI/WIN Polish	CP852 MS-DOS Latin2
Czech	Printer Mode Baud rate: 9600 Bit / Parity: 8 bit None 1 Handshake: Xon/Xoff Language: ANSI/WIN Czech	CP852 MS-DOS Latin2
Hungarian	Printer Mode Baud rate: 9600 Bit / Parity: 8 bit None 1 Handshake: Xon/Xoff Language: ANSI/WIN Hungarian	CP852 MS-DOS Latin2
Portuguese	Printer Mode Baud rate: 9600 Bit / Parity: 8 bit None 1 Handshake: Xon/Xoff Language: ANSI/WIN Portuguese	CP860 MS-DOS Portuguese
English, French, German, Italy, Spanish, Chinese, Japanese, Korean	Printer Mode Baud rate: 9600 Bit / Parity: 8 bit None 1 Handshake: Xon/Xoff	Factory Settings

9 Consumables & Spare Parts

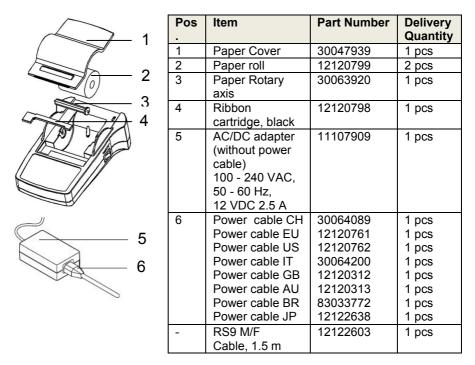

10 RS232C Interface

The SF40A printer is equipped with an RS232C interface to connect OHAUS instruments.

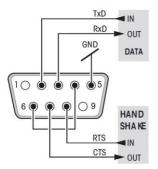

- 9-pin male connector
- Matching to other device (transmission parameters) see chapter 4.

11 Technical Data

Dowor oupply	AC/DC adapter: 12 VDC, 2.5 A		
Power supply	•		
	Primary: 100 – 240 VAC, -50 -60 Hz		
	Voltage fluctuations:		
	up to $\pm 10\%$ of the nominal voltage		
	Secondary: 12 VDC, 2.5 A		
	(with electronic overload protection)		
Cable to AC/DC adapter	3-core, with country-specific plug		
Power line voltage	12 VDC		
Power consumption	Max. 2.5 A		
Altitude	up to 2000 m		
Ambient temperature	0 °C – 40 °C		
Relative humidity	Max. 80% at 31 °C, decreasing linearly to 50% at 40 °C, noncondensing		
Overvoltage category	Class II		
Pollution degree	2		
Printing unit	Dot matrix printer 5x7, 40-character line length		
Printing speed	1.2 lines per second		
Ribbon cartridge	Exchangeable, black		
Paper roll	Standard paper 57.5 mm x ø 50 to 60 mm, integrated in housing, commercial size		
Print quality	Light resistant and thermally stable printing (GLP, GMP, ISO9001)		
Interface	RS232C		
Printer dimensions	(WxDxH) 120 mm x 203 mm x 73 mm		
Packaging dimensions	(WxDxH) 255 mm x 205 mm x 150 mm		
Net weight	740 g (paper roll incl.)		
Compatibility limits	Function Date / Time:		
	not applicable on balances with built-in clock (built-in clock has priority)		
	Functions Zeroing / Taring, Statistics and Totalization:		
	not applicable on all models of balances		



OHAUS Corporation 7 Campus Drive Suite 310 Parsippany, NJ 07054 USA Tel: +1 973 377 9000 Fax: +1 973 944 7177

With offices worldwide / Con oficinas alrededor del mundo / Avec des bureaux dans le monde entier / Weltweite Geshäftsstellen / Con uffici in tutto il mondo.

www.ohaus.com



P/N 30063919 © 2012 OHAUS Corporation, all rights reserved / todos los derechos reservados / tous droits réservés / Alle Rechte vorbehalten / tutti i diritti riservati.

Printed in China / Impreso en la China / Imprimé en Chine / Gedruckt in China / Stampato in Cina